

**<別紙> Cedex Bio  
冷却ユニットの性能劣化に関わる温度制御異常アラーム**

アラームコード

**14.770            Cooling power is too low.**

内容    The cooling power is too low.

処置    Confirm the message and try to continue working.  
          If the message reappears contact Roche Service.

**14.771            Reagent temperature is not being controlled.**

内容    The reagent temperature is not being controlled by the cooling unit.

処置    Confirm the message and try to continue working.  
          If the message reappears, call Roche Service.

**14.772            Fluid cooling pressure is not being controlled.**

内容    The fluid cooling pressure for the cooling unit is not being controlled.

処置    Confirm the message and try to continue working.  
          If the message reappears, call Roche Service.

**14.773            Cooling pressure too low.**

内容    The cooling pressure of the cooling unit is too low.

処置    Confirm the message and try to continue working.  
          If the message reappears, contact Roche Service.

**14.4211           Sensor problems in the cooling unit.**

内容    Sensor problems in the cooling unit.

処置    Confirm the message and try to continue working.  
          If the message reappears, contact Roche Service.